# Electrolux

**MRF 120/35  
MRF 280/35  
MRF 400/35  
MRF 600/35**

**MRF 240/35  
MRF 350/35  
MRF 480/35**

**MRF240/86  
MRF480/86  
MRF401/86  
MRF601/86**

<b>OPERATING INSTRUCTIONS</b>	PAGE 3
<b>BEDIENUNGSANLEITUNG</b>	SEITE 11
<b>NOTICE D' EMPLOI</b>	PAGE 19
<b>GEBRUIKSAANWIJZING</b>	PAGINA 27
<b>ISTRUZIONI PER L' USO</b>	PAGINA 35
<b>INSTRUCCIONES DE MANEJO</b>	PÁGINA 43
<b>MANUAL DE UTILIZAÇÃO</b>	PÁGINA 51
<b>BRUKSANVISNING</b>	SIDA 59
<b>HÅNDBOK</b>	SIDE 67
<b>BETJENINGSVEJLEDNING</b>	SIDE 75
<b>KÄYTTÖOHJE</b>	SIVU 83

# CONTENTS

<b>1</b>	<b>GENERAL WARNINGS .....</b>	<b>4</b>
<b>2</b>	<b>ENVIRONMENTAL PROTECTION .....</b>	<b>4</b>
<b>3</b>	<b>SAFETY .....</b>	<b>4</b>
<b>4</b>	<b>TRANSPORTATION DAMAGES .....</b>	<b>4</b>
<b>5</b>	<b>POSITIONING .....</b>	<b>5</b>
<b>6</b>	<b>CLEANING / DISINFECTION .....</b>	<b>5</b>
<b>7</b>	<b>ELECTRICAL CONNECTION .....</b>	<b>6</b>
<b>8</b>	<b>FUNCTIONING OF THE LOCK .....</b>	<b>6</b>
<b>9</b>	<b>ELECTRONIC SUMMARY .....</b>	<b>6</b>
<b>10</b>	<b>SWITCHING ON .....</b>	<b>6</b>
<b>11</b>	<b>TEMPERATURE DISPLAY .....</b>	<b>7</b>
<b>12</b>	<b>TEMP-SET .....</b>	<b>7</b>
<b>13</b>	<b>ALARM-SET .....</b>	<b>7</b>
<b>14</b>	<b>ALARM-FUNCTION .....</b>	<b>8</b>
<b>15</b>	<b>MEMORY .....</b>	<b>8</b>
<b>16</b>	<b>AC-FAIL .....</b>	<b>8</b>
<b>17</b>	<b>ACCUMULATOR .....</b>	<b>9</b>
<b>18</b>	<b>EXTERNAL ALARM FUNCTION .....</b>	<b>9</b>
<b>19</b>	<b>REASONS FOR ALARM .....</b>	<b>9</b>
<b>20</b>	<b>DEFROSTING .....</b>	<b>9</b>
<b>21</b>	<b>MAINTENANCE .....</b>	<b>10</b>
<b>22</b>	<b>WHEELS .....</b>	<b>10</b>
<b>23</b>	<b>TEMPERATURE RECORDER .....</b>	<b>10</b>

THE EQUIPMENT CONFORMS TO THE EMC DIRECTIVE 89/336/EEC, AS WELL AS TO THE LOW-VOLTAGE DIRECTIVES 73/23/EEC AND 93/68/EEC..

**1****GENERAL WARNINGS**

- Before starting the equipment carefully read the instructions for use, as well as all the work, operation and maintenance safety warnings. Only in this way are correct use and excellent efficiency guaranteed.
- Keep these instructions near the equipment for any further consultation by the various operators regarding the operation and safety standards.
- The equipment must be used exclusively by adults. Keep out of reach of children.
- All the installation and adjustment work on the equipment as well as the maintenance or repair work must be done exclusively by a specialised technician from the Service Centre. The work carried out by people who are not experts can compromise the efficiency of the equipment and cause damage to people or things. Use only original spare parts.
- The modifications made on the equipment can cause damage or injuries.
- Enough space must be allowed around the equipment for air circulation (min. 75 mm) . For this purpose, refer to the installation warnings / instructions.
- The equipment must be put against a wall, usually with the back against the wall, so the parts that are heated are not accessible. Leave enough free space so that good air circulation is guaranteed.
- Make sure that the power supply cable is not stuck or folded during installation or shifting of the equipment.
- Before carrying out the cleaning and maintenance operations disconnect the equipment and remove the plug. Do not pull the wire. If you can not reach the plug disconnect the protection device.
- The cooling system on the rear side and inside the equipment contains a cooler. The equipment and the products preserved in it can undergo remarkable damage if such a system is no longer hermetically sealed. Be careful that blunt objects are not brought in contact with the cooling system.

**2****ENVIRONMENTAL PROTECTION**

- Contribution to the protection of the environment: it is important to appropriately and regularly discharge rubbish.
- The material used for packaging can usually be recycled and must be able to be reused.


**3****SAFETY**

- The products stored are very cold. To protect against burning, gloves should be worn.
- Before scrapping an old refrigerator remove the door so that no children can lock themselves inside while playing.


**4****TRANSPORT**

- The appliance may only be transported while in an upright position (maximum tilt angle 45°).
- Check if the equipment has been damaged. If it is damaged due to transportation IMMEDIATELY turn to the supplier or the dealer showing the delivery note or receipt.
- Equipment that has been damaged by transportation must not be operated!


- The room in which the equipment is to be installed must be dry and suitably aired, avoiding direct exposure to the sun or positioning near a heat source.
- Make sure that the equipment is positioned in a stable manner and on a flat surface. It must never be put next to other equipment. By loosening or tightening the feet, they can be adapted to the desired level.
- For exceptional functioning there must be good airing around the equipment. Make sure that in the lower, rear and top part of the equipment there are no obstacles to air circulation. The free space around the appliance must be at least 75 mm.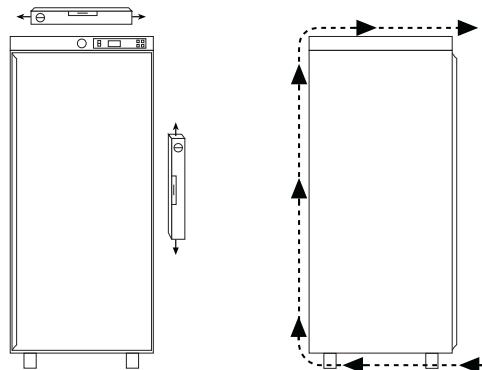


- Models MRF400/35 and MRF600/35 should be attached to the wall using the catch provided, to ensure they do not overturn.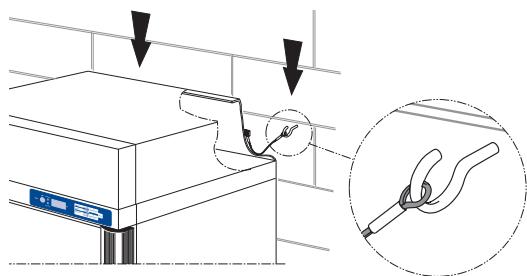

- Models MRF400/35 and MRF600/35 are fitted with an automatic defroster. To ensure optimum defrosting, the built-in defrosting dial should be set to the present time before commissioning.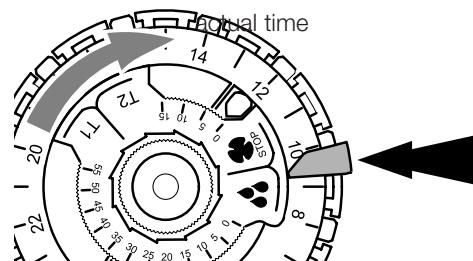


- The door handle on models MRF401/86 and 601/86 has been moved by 90° at the factory end to facilitate packaging and transporting, and should be turned during installation in accordance with the diagram across.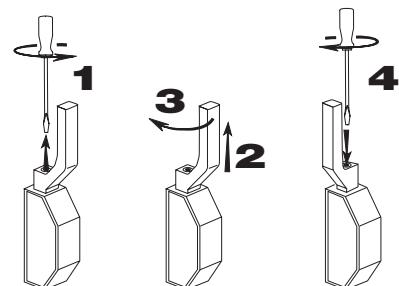

- Before using clean or disinfect all the internal and external parts. Use only delicate detergents. Completely avoid using aggressive, corrosive or abrasive detergents, steel wool, abrasive sponges or chemical solvents. Be careful that during cleaning, liquids do not filter into the ventilator housing.
- If the internal equipment is made of steel, there are plastic foils on different parts of the doors and drawers. Remove these foils before starting the refrigerator.

- Before carrying out the electrical connection of the equipment, make sure that the data stated on the data plate (in the top left on the internal container or on the back side) correspond to the local values. It is important to connect the equipment to a ground clamp.
- Before connecting the equipment and activating it for the first time, it must be in a vertical position for about 30 minutes. If the equipment was transported horizontally, wait 4 hours.
- To avoid the equipment not functioning due to any breakdowns to other electrical equipment it is necessary to be connected to a separate electrical circuit. In no case must it be connected together with other electrical equipment by means of a multiple plug, to a single electrical outlet.

- When the key is in the horizontal position the door is closed and the key can be removed.
- If the unit is to be left permanently open and accessible to all (daytime operation), turn the key 90° to the right to the vertical position and remove it.
- Do not store the keys in reach of children or near the refrigerator.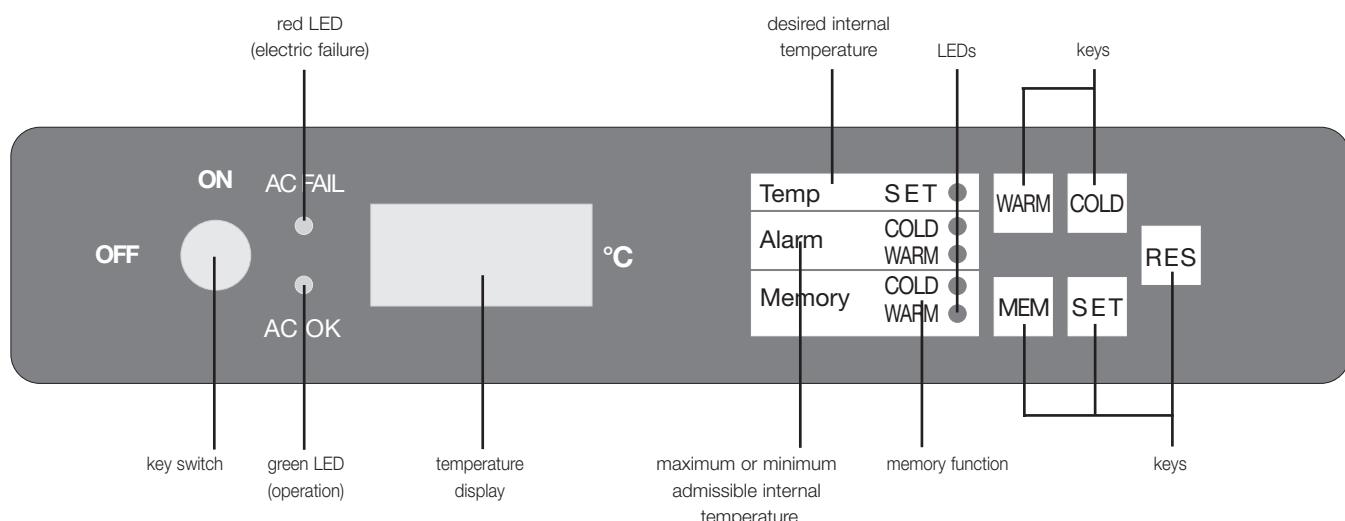

- To operate the equipment turn the key switch on the command panel to the „ON“ position. In this position the key can be taken out to avoid the equipment from disconnecting by mistake.
- Upon ignition all the LEDs light up constantly for about 5 seconds. The acoustic alarm is simultaneously activated and the following temperature value appears on the screen **-18,8**.
- Following this initialization only the green LED remains lit. It indicates that the equipment is connected and that it is live. The present temperature inside the refrigerator appears on the screen now and corresponds to room temperature after ignition. As the display shows temperatures in a particular range only, temperatures below this range are shown with **--** and above with **---**.
- After the first ignition check the correct functioning of the equipment and if the set temperature is reached after 24 hours.
- Upon the first ignition the alarm remains inactive until the internal temperature reaches the set value. Only after this occurs is the alarm active.
- When switching the appliance off with the key switch, all parameters of the electronics are set back to factory settings.

## 11

## TEMPERATURE DISPLAY


- On the display screen the internal refrigerator temperature is displayed.

- For /35 appliances:**

If the temperature is colder than -60°C, the display  will appear; temperatures warmer than -10°C will be shown with .

- For /86 appliances:**


If the temperature is colder than -99°C, the display  will appear; temperatures warmer than -60°C will be shown with .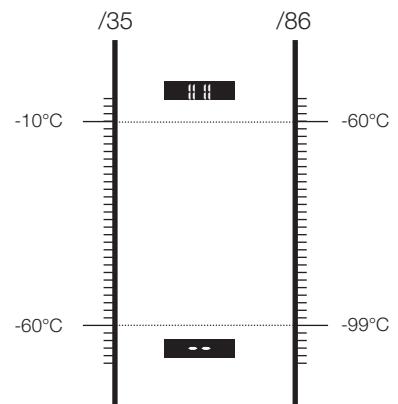

## 12

## TEMP-SET



- By „Temp-Set“ we mean the desired internal temperature of the refrigerator. This temperature has been preset in the factory. If needed it can be adjusted. Such adjustment must be done by a technician from the Service Centre. If the „Temp-Set“ is changed, the „Alarm-Sets“ change automatically.
- In order to be able to detect the „Temp-Set“ set, press the „SET“ key. The „Temp-Set“ LED will light up and the „Temp-Set“ value set will be displayed as long as the „SET“ key remains pressed.
- Pre-set values:
  - For /35 appliances:** -35°C
  - For /86 appliances:** -82°C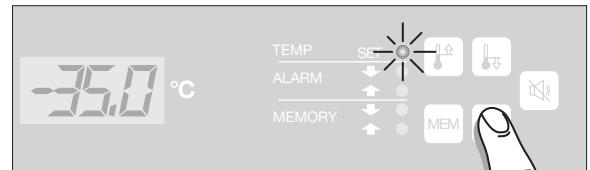


## 13

## ALARM-SET



- By „Alarm-Set“ we mean the maximum and minimum admissible internal temperature. It can be modified as follows, according to needs :
  - For /35-models :** Alarm COLD = minimum value set: -55°C , maximum value set: 5° below the „Temp-Set“. Alarm WARM = maximum value set: -15°C, minimum value set: 5° above the „Temp-Set“.
  - For /86-models :** Alarm COLD = minimum value set: -95°C, maximum value set: 5° below the „Temp-Set“. Alarm WARM = maximum value set: -65°C, minimum value set: 5° above the „Temp-Set“.
- By pressing the relative WARM and COLD keys the „Alarm-Set“ values set can be detected. The relative „WARM-alarm“ or „COLD-alarm“ LEDs light up while the set temperature is displayed. The following values have been preset:
  - For /35-models :** Alarm COLD : -40°C      Alarm WARM : -30°C
  - For /86-models :** Alarm COLD : -87°C      Alarm WARM : -77°C
- In the setting of the „Alarm-Set“ proceed as follows:
  - Alarm-Cold = Press the “SET” and “COLD” key at the same time; the “Alarm-Cold” LED lights up and the temperature changes in 1°C stages from the set value up to the end value. Once this value has been reached, the display moves back to the original value. In order to set a value, simply release the keys once the temperature required is displayed.
  - Alarm-Warm = Press the “SET” and “WARM” key at the same time; the “Alarm-Warm” LED lights up and the temperature changes in 1°C stages from the set value up to the end value. Once this value has been reached, the display moves back to the original value. In order to set a value, simply release the keys once the temperature required is displayed.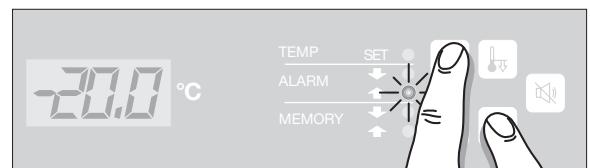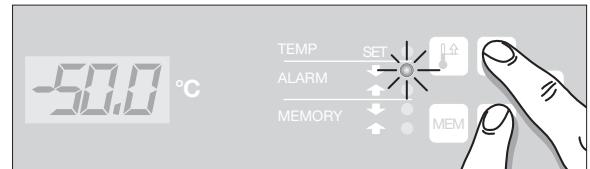


- The measurement of the internal refrigerator temperature is carried out by means of 2 sensors placed inside the refrigerator. The temperature indicated comes from the average value of the two temperatures measured. To prevent the air temperature being measured in this case instead of the temperature of the goods being stored, the two sensors are embedded in a reference unit which is filled with liquid. This simulates the temperature time lag found in the goods.
- The temperature alarm is activated when :
  - the average value of the two temperatures detected does not fall in the set area (COLD alarm - WARM alarm),
  - one of the two detected temperatures does not fall in the set area even if the average value is within the acceptable margin,
  - both of the values do not fall in the margin
- When the alarm starts there is an audible alarm and the respective alarm-LED starts to flash (Alarm-COLD or Alarm-WARM). On the display the temperatures of the two sensors appear successively at regular second intervals.

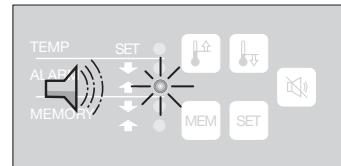
- By pressing the „RES“ key the audible alarm is discontinued, the flickering light display of the two temperature values indicated by the sensors continues as well as the flashing function of the alarm LED until the temperature irregularity is eliminated. If this does not happen within 20 minutes the audible alarm goes off again.
- The alarm turns off automatically as soon as the temperature goes back to the normal values. To warn the operator of the fact that the alarm has gone off, the LED-Mem continues to flash until the temperature is read by means of the Memory function.

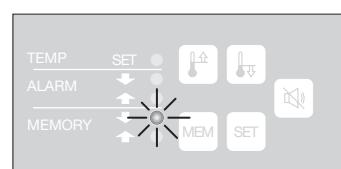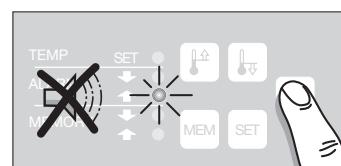- The maximum and minimum temperature shown after the last ignition can be determined with the help of the „Memory“ function. For this purpose press the „MEM“ and „WARM“ keys simultaneously for the maximum temperature and „MEM“ and „COLD“ for the minimum temperature. The „Memory-COLD“ or the“Memory-Max“ LED turns on displaying the relative value.
- To cancel the memorized values press the „MEM“ and „RES“ keys simultaneously. The „Memory“ LEDs turn on and the two values are cancelled.

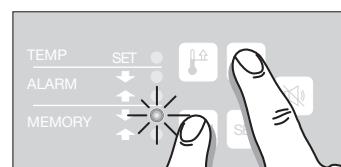- If the voltage should be discontinued (for having disconnected the plug or due to a breakdown in the safety valve), the green LED turns off on the command panel and the red LED begins to blink. Furthermore, the audible alarm is activated. It can be deactivated temporarily with the „RES“ key.
- During the „AC-fail“ phase no temperature is displayed. The temperature measurement continues and the maximum and minimum values can be recalled as soon as the network voltage is present (see „Memory“).
- When the network voltage is missing the set data cannot be modified or cancelled.
- If during this phase there is a temperature increase the „WARM-Alarm“ LED as well as the red „AC-FAIL“ LED will flash.
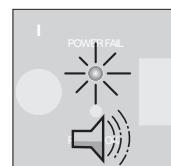
## 17


## ACCUMULATOR

- In the electronic part an accumulator is included. If there is a lack of electricity, it carries out the alarm and temperature measurement functions for a maximum of 22 hours.
- This accumulator is charged automatically, as soon as the equipment plug is connected.
- The accumulator should be replaced every three years. The replacement must be carried out only by the Technical Service Centre.

## 18

## EXTERNAL ALARM FUNCTION

- On the back of the electronic circuit there are two potential-free outlets that can be used to command an external alarm system. These outlets are guided to two sockets at the rear of the refrigerator where an external optical or acoustical alarm can be connected.
- The external alarm function only reports a temperature alarm. **An AC-Fail alarm is not indicated.**
- The two potential-free outlets are related to a make-contact.
- The connected voltage must not exceed 30V DC. The current is limited to 2A.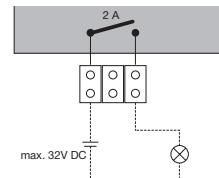

## 19

## ALARM REASONS


- As already previously described two types of alarm can be distinguished: a temperature alarm and the „AC-Fail“ alarm.
- The sounding of the AC-fail alarm can happen for the following reasons:
  - plug disconnected from the outlet
  - outlet without electricity (faulty safety valve, lack of electricity, etc.)
- When there is a temperature alarm, check for the following possible causes:
  - Has the door been left open longer than necessary, causing an increase in the internal temperature?
  - Have too many hot or cold products been placed in the refrigerator at the same time?
  - Have products with packages that have been heated been placed in the refrigerator?
  - Has a layer of ice formed on the evaporator plate (at the back inside)? This would mean that the defrosting does not work and that the equipment is not able to supply sufficient cooling power.
  - Is the room temperature too high, that is, over 32°C?
- If the cause of the alarm cannot be immediately eliminated, there is the danger that the products become too hot or too cold. In this case it is recommended to empty the refrigerator and put the contents in other equipment.
- Check the functioning of the refrigerator after having eliminated all the described causes. Warn the Assistance Service if after a certain period of time the temperature values do not go back to normal.

## 20

## DEFROSTING

- Excessive frosting-over of the evaporator prevents the appliance's cold chamber from refrigerating and reduces refrigeration capacity. Regular defrosting is therefore essential.
- Models MRF400/35 and MRF600/35 are fitted with an automatic defroster. To ensure optimum defrosting, the built-in defrosting dial should be set to the present time before commissioning and after the appliance has been switched off for a long time (over one hour).
- All other appliances should be defrosted manually once a layer of ice over 1 cm thick has formed.
- For defrosting, proceed as follows:  
Re-store products in another appliance. Set the locking switch to the 'OFF' position and, with the door open, wait until the ice has melted.
- After defrosting, the inside of the appliance should be thoroughly cleaned and dried.


- To clean the refrigerator use delicate detergents. Completely avoid the use of aggressive or corrosive products, abrasive powder, steel wool, abrasive sponges or chemical solvents.
- The condenser (if not lathered) should be dusted twice per year with a hand brush or a vacuum cleaner. The condenser is located at the rear end or under the cover, depending on the model.
- For /86 appliances, a filter is located behind the ventilation grille, which should be regularly dusted with a hand brush or a vacuum cleaner.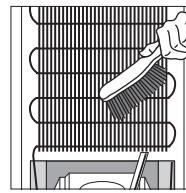


- For MRF400/35 and MRF600/35 :**

As an option, these appliances may come provided with 4 mobile transport rollers. In this case, both of the front rollers will be fitted with a locking device.

- To lock the brake push the pedal towards the bottom and to unlock it push the smallest lever forward.
- Warning :** The brake must be locked or unlocked using only feet.
- Appliances MRF240, MRF480, MRF401 and MRF601 may be safeguarded from moving away by adjusting both locking screws.

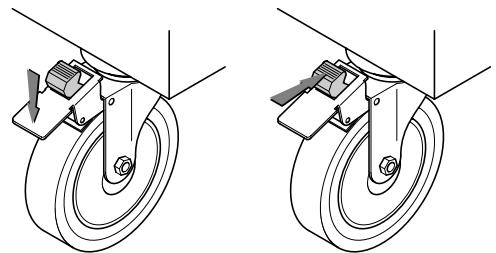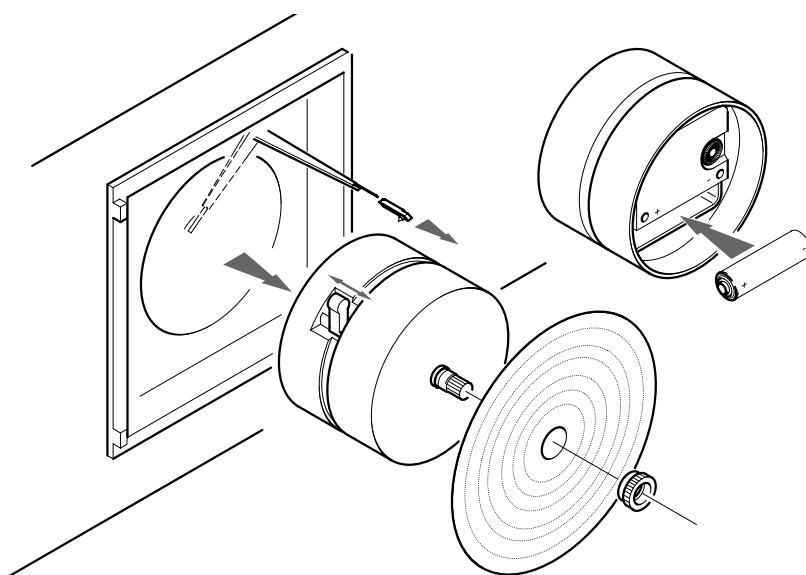- By means of the temperature recorder for a certain time interval the course of the temperature is recorded. The time can be programmed according to needs, at 24 hours or 7 days.
- Before starting up, it is recommended to check the set time in the following manner: raise the lever with the pen point, loosen the lock nut on the recorder axis and remove the recorder disk. Pull the recorder axis and remove the attachment from the housing. On the side of the attachment there is the switch with which you can set the desired time.
- The disks for the two time intervals are different from each other. Check if the suitable disk has been inserted.
- The recorder horology is fed by a 1.5 V battery which is found in the lower part of the recorder attachment.
- To replace the pen point, remove the old one from the bracket and insert the new one until it stops. Avoid touching the point with your hands.

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